

PRODUCT BULLETIN: Cellometer Auto

Cellometer Auto Compatibility with Macintosh.



Cellometer Auto can be easily run on a Mac meeting the following minimum requirements:

- 2.0 GHz or higher Intel based Mac running OS X Leopard, with 1 GB RAM.
- Parallels Desktop 3.0 running Windows XP (www.parallels.com)



After launching Parallels, simply load Cellometer Software and follow standard installation instructions. The Cellometer USB 2.0 cable must be connected after Parallels is launched to ensure the instrument is registered in the Windows environment.

Consult Parallels documentation for more information about transferring output images and data files to the Mac environment.

For more information, please contact us at: support@nexcelom.com, or call 1-978-327-5340.